




SIGMOIDA ANALYSIS SOFTWARE

- Automated result interpretation
- Reporting in 2 seconds
- The **offline** software ensures data safety
- Automated report transfer to the Laboratory Information Systems.
- Human error elimination
- Ready-to-use for multiple kits

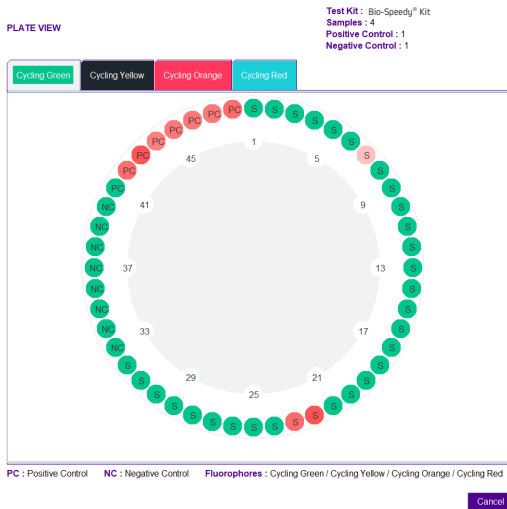

Mic qPCR Cycler Plate Setup Interface

PLATE VIEW

Test Kit : Bio-Speedy® Kit
Samples : 40
Positive Control : 4
Negative Control : 4

| | FAM | HEX | ROX | Cy5 | 1 | 2 | 3 | 4 | 5 | 6 | 7 | 8 | 9 | 10 | 11 | 12 |
|---|-----|-----|-----|-----|---|---|---|---|---|---|---|---|---|----|----|----|
| A | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| B | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| C | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| D | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| E | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| F | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| G | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |
| H | | | | | S | S | S | S | S | S | S | S | S | S | NC | PC |

PC : Positive Control NC : Negative Control Fluorophores : FAM / HEX / ROX / Cy5

Bio-Rad Plate Setup Interface

Report

Test Kit : Bio-Speedy® Kit
File : CVD FLU DATA 1.pcrd
Date : 03 Feb 2023 09:10:22
User : Admin
Version : V 2.1.3

Show in plate view

All (171) Positive (37) Negative (101) Issues (33) Controls (39)

Select All

| | Well | Sample | Target | Cq | Result | Warnings |
|-------------------------------------|------|---------------|-------------|-------|----------|----------|
| <input checked="" type="checkbox"/> | A1 | CVD HASTA 1 | SARS-CoV-2 | 22.32 | Positive | |
| <input checked="" type="checkbox"/> | A1 | CVD HASTA 1 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | A1 | CVD HASTA 1 | Influenza A | N/A | Negative | |
| <input checked="" type="checkbox"/> | B1 | CVD HASTA 1 | SARS-CoV-2 | 24.04 | Positive | |
| <input checked="" type="checkbox"/> | B1 | CVD HASTA 1 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | B1 | CVD HASTA 1 | Influenza A | N/A | Negative | |
| <input checked="" type="checkbox"/> | D1 | INF A HASTA 1 | SARS-CoV-2 | N/A | Negative | |
| <input checked="" type="checkbox"/> | D1 | INF A HASTA 1 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | D1 | INF A HASTA 1 | Influenza A | N/A | Negative | |
| <input checked="" type="checkbox"/> | E1 | INF B HASTA 1 | SARS-CoV-2 | N/A | Negative | |
| <input checked="" type="checkbox"/> | E1 | INF B HASTA 1 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | E1 | INF B HASTA 1 | Influenza A | N/A | Negative | |
| <input checked="" type="checkbox"/> | A2 | CVD HASTA 2 | SARS-CoV-2 | 25.2 | Positive | |
| <input checked="" type="checkbox"/> | A2 | CVD HASTA 2 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | A2 | CVD HASTA 2 | Influenza A | N/A | Negative | |
| <input checked="" type="checkbox"/> | B2 | CVD HASTA 2 | SARS-CoV-2 | 24.81 | Positive | |
| <input checked="" type="checkbox"/> | B2 | CVD HASTA 2 | Influenza B | N/A | Negative | |
| <input checked="" type="checkbox"/> | B2 | CVD HASTA 2 | Influenza A | N/A | Negative | |

Result Reporting Table



Head Office / Manufacture Office : Huzur Mah. Metin Oktay Cad. Nurol Life Sitesi D Blok No:3/31 Sarıyer-Istanbul-TURKEY
R&D / Project Office : Büyükdere Cad. No: 237 Nörâmin İş Merkezi Kat: B4 Maslak Sarıyer-Istanbul-TURKEY
 Phone: +90 212 285 10 17 - www.bioeksen.com - e-mail: info@bioeksen.com.tr